

# 7-Wire CerePlex to Plastics1 Commutator Cable

*PN 7910*



## **Blackrock Microsystems, LLC**

630 Komas Drive, Suite 200

Salt Lake City, UT 84108


Tel: (866) 806-3692

[www.blackrockmicro.com](http://www.blackrockmicro.com)

# 7-Wire CerePlex to Plastics1 Commutator Cable - Instructions for Use

## 1.1 7-Wire CerePlex to Plastics1 Commutator Cable

The 7-Wire CerePlex to Plastics1 Commutator Cable provides an interface between the outputs of a Blackrock CerePlex M32 Headstage and the bottom of an 18-Channel Plastics1 Commutator. The pin-out for this cable is provided in the figures below. Pin-outs are viewed looking into the connector.

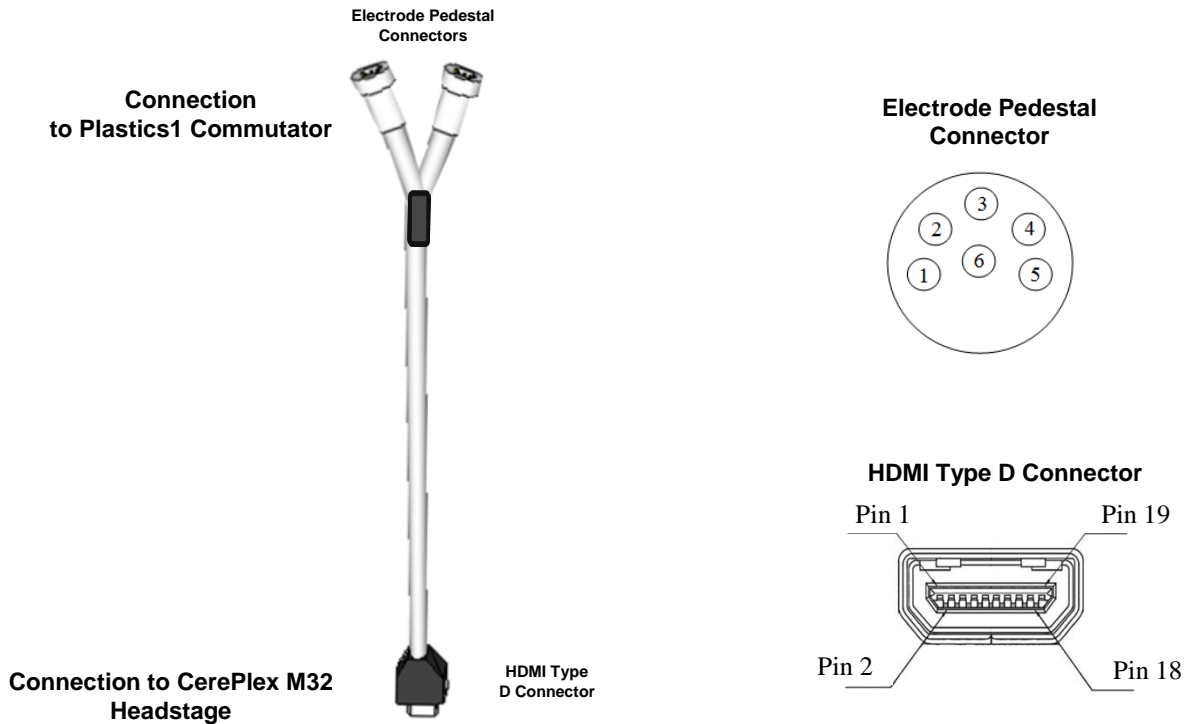

**Electrode Pedestal Connector #1 (WHITE)**

Electrode Pedestal Connector Pin	HDMI Type D Connector Pin
1	1 (Ground)
2	19 (V+)
3	16 (V-)
4	Not Connected
5	Not Connected
6	Not Connected

**Electrode Pedestal Connector #2 (BLACK)**

Electrode Pedestal Connector Pin	HDMI Type D Connector Pin
1	9 (DATA+)
2	11 (DATA-)
3	12 (+CLK)
4	14 (-CLK)
5	Not Connected
6	Not Connected

**\*\*NOTE:** Connectors are viewed looking into the connector.

## 2 Warranty

Blackrock Microsystems, LLC warrants that its products are free from defects in materials and manufacturing for a period of one year from the date of shipment. Blackrock will, at its option, repair or replace any product that does not comply with this warranty. This warranty is voided by:

1. Any modification or attempted modification to the product done by anyone other than an authorized Blackrock employee
2. Any abuse, negligent handling or misapplication of the product.

This constitutes the sole warranty made by Blackrock, LLC. There are no other warranties, expressed or implied, which extend beyond those described herein or to anyone other than the original purchaser, including the implied warranties of merchantability and fitness for a particular purpose. In no event shall Blackrock Microsystems, Inc. be liable for any incidental or consequential damages, or for the infringement of any patent rights or third party rights due to the use of its products.

### 2.1 Return Merchandise Authorization (RMA)

In the unlikely event that your adaptor needs to be returned to Blackrock for service or maintenance, do not send any equipment back without a Return Merchandise Authorization Number. An RMA number will be issued to you by a Blackrock representative. If you need to obtain an RMA number, you may contact a product support representative at (801) 582-5533 or toll free at (866) 806-3692.

Once an RMA number has been issued, it is important to safely pack the returned item for shipping back to Blackrock. It is preferred that you save the original boxes and packing materials that your system arrived in for return shipment. Please address the package as follows:

**Blackrock Microsystems, LLC**  
**ATTN: RMA#**  
**630 Komas Drive, Suite 200**  
**Salt Lake City, UT 84108 USA**  
**Tel: (801) 582-5533**