


Blackrock QHA-32 / QHA-2x16 Adapter

PN-8808 & PN-8809

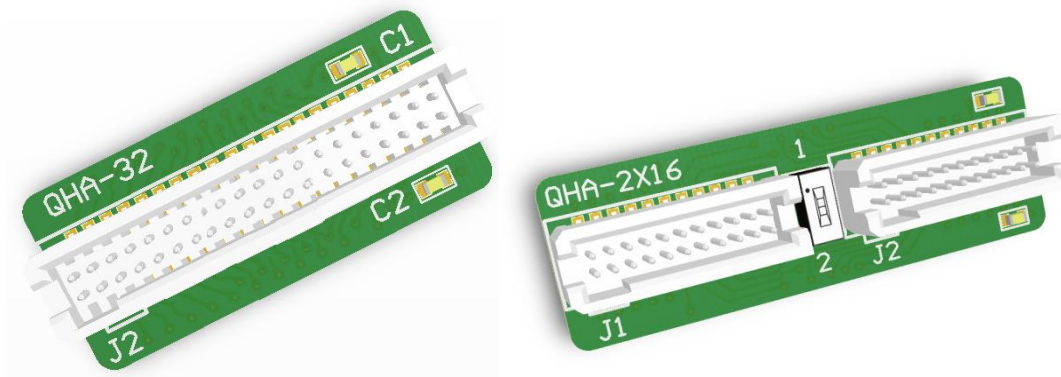

Blackrock Microsystems, LLC

630 Komas Drive • Suite 200

Salt Lake City • UT 84108

T: +1 (801) 582-5533

www.blackrockmicro.com

support@blackrockmicro.com

Electrophysiology, Simplified.

1 Introduction

The QHA adapter provides an interface for connecting Blackrock's legacy connector to the new CerePlex A.

2 Hardware

2.1 QHA-32 Adapter Pin-out

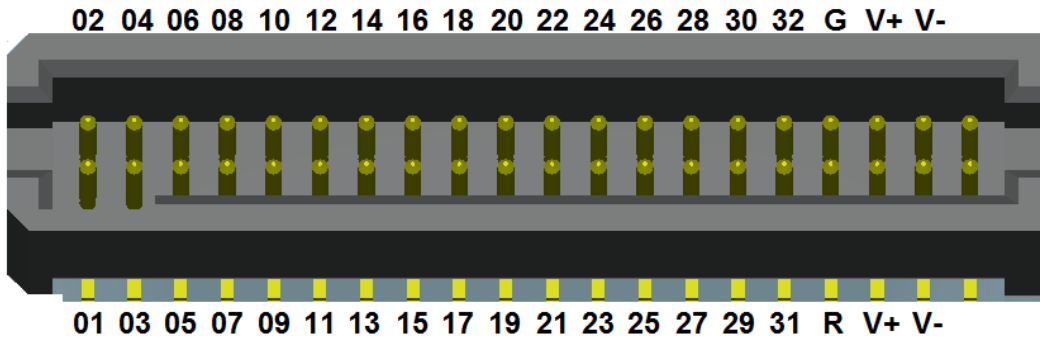


Figure 2-1: Blackrock QHA-32 Adapter Pin-Out

2.2 QHA-2x16 Adapter Pin-out


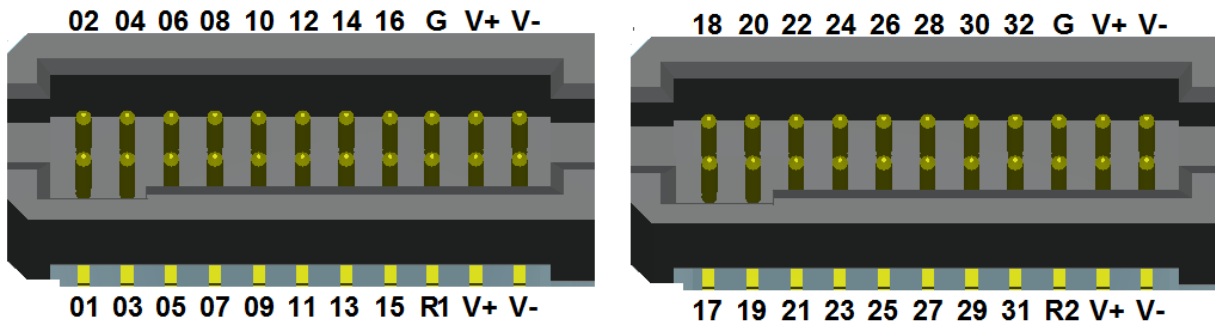


Figure 2-1: Blackrock QHA-2x16 Adapter Pin-out

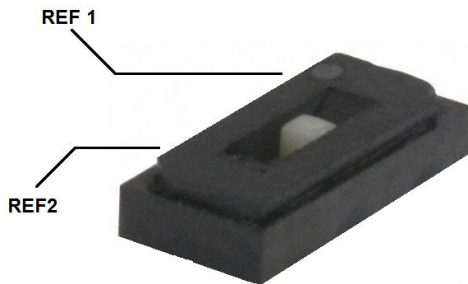

The switch on the QHA-2x16 is for reference selection. If the switch is in the top position, both of the connectors will use the reference R1 from the left connector. If the switch is in the bottom position, both of the connectors will use the reference R2 from the connector on the right.

Figure 2-3: Blackrock QHA-2x16 Adapter reference selection.

3 Warranty

Blackrock Microsystems warrants that its products are free from defects in materials and manufacturing for a period of one year from the date of shipment. Blackrock will, at its option, repair or replace any product that does not comply with this warranty. This warranty is voided by:

1. Any modification or attempted modification to the product done by anyone other than an authorized Blackrock employee
2. Any abuse, negligent handling or misapplication of the product.

This constitutes the sole warranty made by Blackrock, There are no other warranties, expressed or implied, which extend beyond those described herein or to anyone other than the original purchaser, including the implied warranties of merchantability and fitness for a particular purpose. In no event shall Blackrock Microsystems, be liable for any incidental or consequential damages, or for the infringement of any patent rights or third party rights due to the use of its products.

3.1 Return Merchandise Authorization (RMA)

In the unlikely event that your adapter needs to be returned to Blackrock for repair or maintenance, do not send any equipment back without a Return Merchandise Authorization Number. An RMA number will be issued to you by a Blackrock representative. If you need to obtain an RMA number, you may contact a product support representative at (801) 582-5533 or toll free at (866) 806-3692.

Once an RMA number has been issued, it is important to safely pack the returned item for shipping back to Blackrock. It is preferred that you save the original boxes and packing materials that your system arrived in for return shipment. Please address the package as follows:

Blackrock Microsystems
ATTN: RMA#
630 Komas Drive, Suite 200
Salt Lake City, UT 84108 USA
Tel: +1 (801) 582-5533